

What's new... Tainlab WinTec Version 5

Tainlab

Simple & versatile data-gathering systems

New table format for text results...

Text results are now displayed in an improved spreadsheet style table, allowing you to tailor the presentation of your results.

It's easy to hide columns so that you can focus on the data of most interest. Just click the right mouse button and select **Hide**

	Time	Water	Sodium Thiosulphate
	minutes	(°C)	(°C)
1	0.00	20.7	21.5
2	1.00	21.1	21.4
3	2.00	20.7	21.1
4			
5			
6			
7			
8			
9			
10			
11			

New flexibility in Export and Printing...

Select portions of the data, the frequency (eg. every 2nd, every 5th reading), show row numbers, include experiment information and show hidden columns. You can even preview the contents before saving or printing.

Data export made easy...

Just click the right mouse button and select...

- or **Copy all results to clipboard** then paste data directly into Microsoft Excel.
- or use the mouse to select the data you are most interested in, click the right mouse button and select **Copy selected results to clipboard**
- or **Export** to save your results to any of these file formats...
.XLS, .DOC, .CSV, .HTML, .TXT (Tab delimited), Casio calculator .CAT files, Vernier Graphical Analysis or Texas calculator file (CSV - Matrix)


New Printing Facilities...

The new print menu contains all printing options on a single page, so it is easy to format your results tables and graph.

Click the **Print Preview** tab to see how your results will look.

Print Options

- Select Graph or Text results.
- Select Printer and page orientation.
- Add Headers and Footers.
- Show or Hide Table Columns.
- Change Font Size/Text Alignment
- Select a Range to Print.
- Adjust graph size (scaling) and page margins

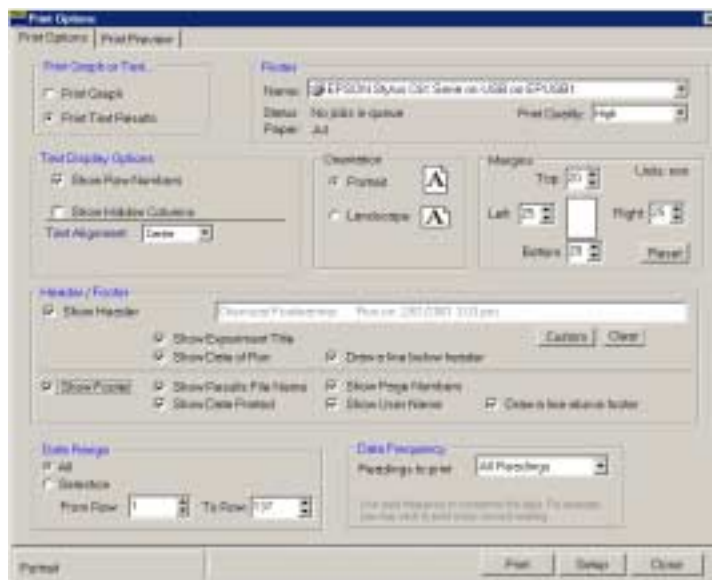

Tainlab software updates are now available via the web

Software Updates are now available for WinTec, ECG, Colourimeter, Tainlab Psychology, Primary Science, PC Digital CRO and Driver Reaction Tester.

Just go to www.tainlab.com and click on the **Software Updates** link.



Tain Electronics Pty Ltd
10 Rowern Court, Box Hill North,
Victoria 3129, Australia
ABN: 70 006 496 706

Tel: 03 9898 7366
Fax: 03 9899 7992
web: www.tainlab.com
email: info@tain.com.au